


Residual Solvents

Column: ZB-624
Dimensions: 60m x 0.32mm x 1.80µm
Flow: Constant flow of 1.0 mL/min of He
Injector: 220°C; 0.6 µL at 20:1 split
Oven: 40°C for 8 min. to 90°C at 8°C/min. to 210°C at 15°C for 8 min.
Detector: MS scanning from 15 to 450 amu
Analytes: All analytes were 140ng on column except

Common Solvents on ZB-624, 60m x 0.32mm x 1.80µm

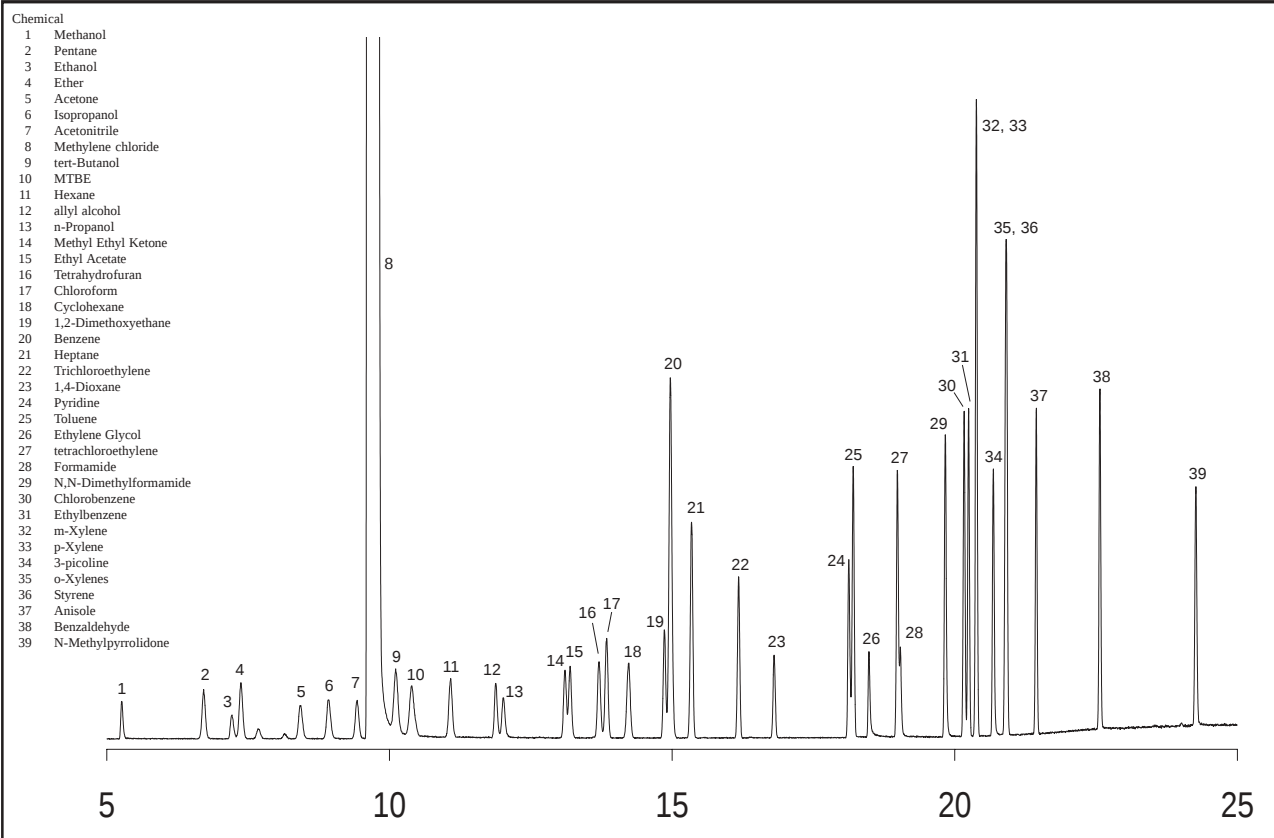

Ordering Information

Order Number	Description
7KM-G005-31	ZB-624 - 60m x 0.32mm x 1.80µm



www.phenomenex.com

Phenomenex products are available worldwide. For the distributor in your country, contact Phenomenex USA, International Department by telephone, fax or e-mail: international@phenomenex.com.

phenomenex
...breaking with tradition

USA	Puerto Rico	Canada	United Kingdom	Germany	New Zealand	Australia
tel.: (310) 212-0555	(800) 541-HPLC	(800) 543-3681	01625-501367	06021-58830-0	09-4780951	1800-553-929
fax: (310) 328-7768	(310) 328-7768	(310) 328-7768	01625-501796	06021-58830-11	09-4780952	1800-553-923
email: info@phenomenex.com	info@phenomenex.com	info@phenomenex.com	ukinfo@phenomenex.com	anfrage@phenomenex.com	info@phenomenex.co.nz	info@phenomenex.com.au