


Nitrogen and Phosphorous Pesticides - EPA Method 507

Column:	Zebtron ZB-5ms	Gold Seal:	Phenomenex single groove (AG0-7518)
Dimensions:	30meter x 0.25mm x 0.25µm	Liner:	Split/Splitless Single Taper (AG0-7516)
Order No.:	7HG-G010-11	Septum:	Phenomenex PhenoRed-400 (AG0-4690)
Injection:	Split 20:1 @ 250°C, 1µL		
Carrier Gas:	Constant Flow Helium 1.0mL/min		
Oven Program:	100°C to 170°C at 20°C/min to 210°C at 4°C/min to 300°C at 20°C/min for 4 min		
Detector:	MSD; 30-400amu		

Notes: Analytes are 50 ppm. Peak identifications are listed on Page 2.

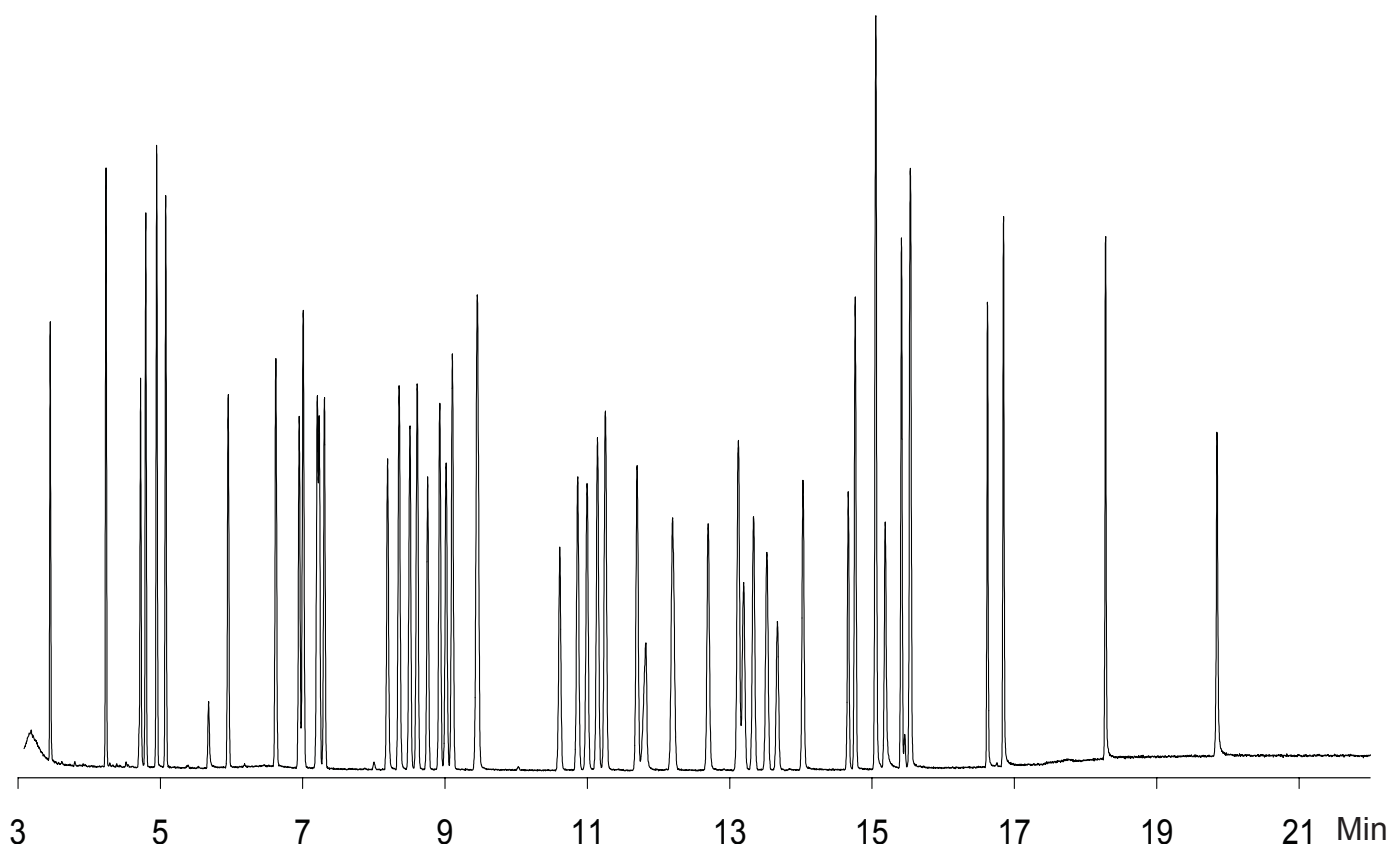


Phenomenex products are available worldwide. For the distributor in your country, contact Phenomenex USA, International Department by telephone, fax or e-mail: international@phenomenex.com.

phenomenex
...breaking with tradition™


USA	Puerto Rico	Canada	United Kingdom	Germany	New Zealand	Australia
mail: 411 Madrid Avenue Torrance, CA 90501-1430 USA	273 Sierra Morena, Suite #104 San Juan, Puerto Rico 00926	411 Madrid Avenue Torrance, CA 90501-1430 USA	Queens Avenue, Hurdfield Ind. Est., Macclesfield, Cheshire SK10 2BN, England	Zeppelinstr. 5 63741 Aschaffenburg Deutschland	P.O. Box 31-601 Milford Auckland New Zealand	P.O. Box 4084 Lane Cove, NSW 1595 Australia
tel.: (310) 212-0555 fax: (310) 328-7768 email: info@phenomenex.com	(800) 541-HPLC (310) 328-7768 info@phenomenex.com	(800) 543-3681 (310) 328-7768 info@phenomenex.com	01625-501367 01625-501796 ukinfo@phenomenex.com	06021-58830-0 06021-58830-11 anfrage@phenomenex.com	09-4780951 09-4780952 info@phenomenex.co.nz	02-9428-6444 02-9428-6445 info@phenomenex.com.au

Nitrogen and Phosphorous Pesticides - EPA Method 507 (continued)

Sample:	1. Dichlorvos	9. Propachlor	17. Atrazine
	2. EPTC	10. Ethoprop	18. Propazine
	3. Mevinphos	11. Cycloate	19. Profluralin
	4. Butylate	12. Chlorpropham	20. Terbufos
	5. Vernolate	13. Trifluralin	21. Pronamide (Propyzamide)
	6. Pebulate (Tillam)	14. Benefin (Benfluralin)	22. Diazinon (Dimpylate)
	7. Tebuthiuron	15. Atraton	23. Disulfoton
	8. Molinate	16. Prometon	24. Terbacil

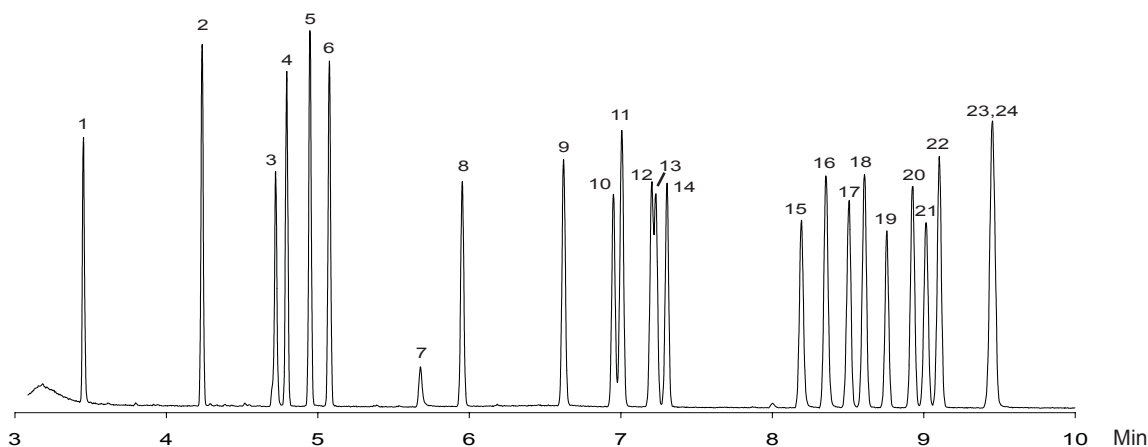

Sample:	25. Metribuzin	34. Diphenamid	43. Fenamiphos
	26. Alachlor	35. MGK-264	44. Tricyclazole
	27. Simetryn	36. Isopropalin	45. Oxadiazon
	28. Ametryn	37. Pendimethalin	46. Oxyfluorfen
	29. Prometryn	38. MGK-264 (Isomer)	47. Carboxin
	30. Terbutryn	39. Merphos	48. Norflurazon
	31. Bromacil	40. Stirofos (Tetrachlorovinphos)	49. Hexazinone
	32. Metolachlor	41. Butachlor	50. Fenarimol
	33. Triadimefon	42. Napropamide	51. Fluridone

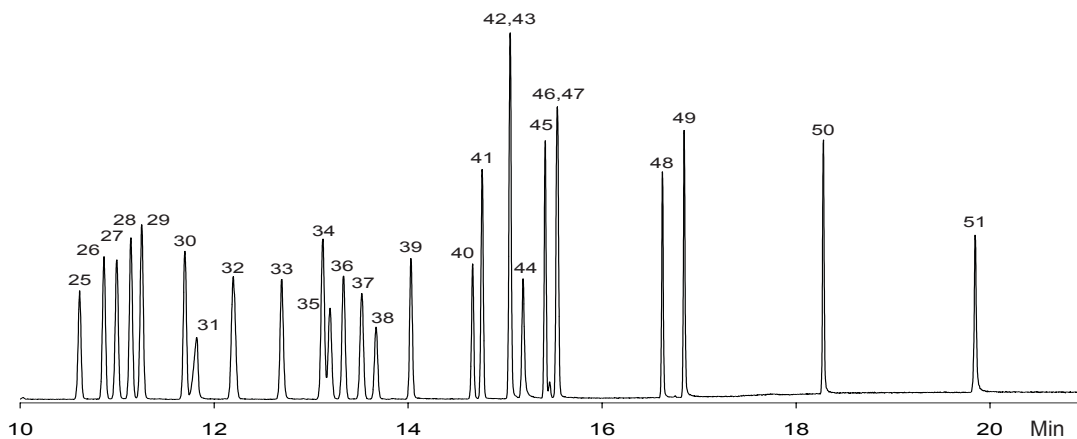

Phenomenex products are available worldwide. For the distributor in your country, contact Phenomenex USA, International Department by telephone, fax or e-mail: international@phenomenex.com.

phenomenex®
...breaking with tradition™

USA mail: 411 Madrid Avenue Torrance, CA 90501-1430 USA tel.: (310) 212-0555 fax: (310) 328-7768 email: info@phenomenex.com	Puerto Rico 273 Sierra Morena, Suite #104 San Juan, Puerto Rico 00926 (800) 541-HPLC (310) 328-7768 info@phenomenex.com	Canada 411 Madrid Avenue Torrance, CA 90501-1430 USA (800) 543-3681 (310) 328-7768 info@phenomenex.com	United Kingdom Queens Avenue, Hurdfield Ind. Est., Macclesfield, Cheshire SK10 2BN, England 01625-501367 01625-501796 ukinfo@phenomenex.com	Germany Zeppelinstr. 5 63741 Aschaffenburg Deutschland 06021-58830-0 06021-58830-11 anfrage@phenomenex.com	New Zealand P.O. Box 31-601 Milford Auckland New Zealand 09-4780951 09-4780952 info@phenomenex.co.nz	Australia P.O. Box 4084 Lane Cove, NSW 1595 Australia 02-9428-6444 02-9428-6445 info@phenomenex.com.au
--	--	---	--	---	---	---