


USP 467 - Method V

App ID 14799

Column: Zebtron ZB-624
Dimensions: 30meter x 0.53mm x 3.0µm with 5m Z-Guard
Order No.: 7HK-G005-36; 7AK-G000-00-GZ0
Injection: Split 2:1 @ 140°C, 1uL
Carrier Gas: Constant Flow Helium 5.0 mL/min
Oven Program: 35°C for 20 min to 240°C at 35°C/min
Detector: FID @ 260°C
Sample:

1. Methylene Chloride
2. Chloroform
3. Benzene
4. Trichloroethylene
5. 1,4-Dioxane

Notes: Methylene Chloride-10ug/mL; Chloroform-1ug/mL; Benzene, 1,4-Dioxane, and Trichloroethylene at 2 ug/mL in water.

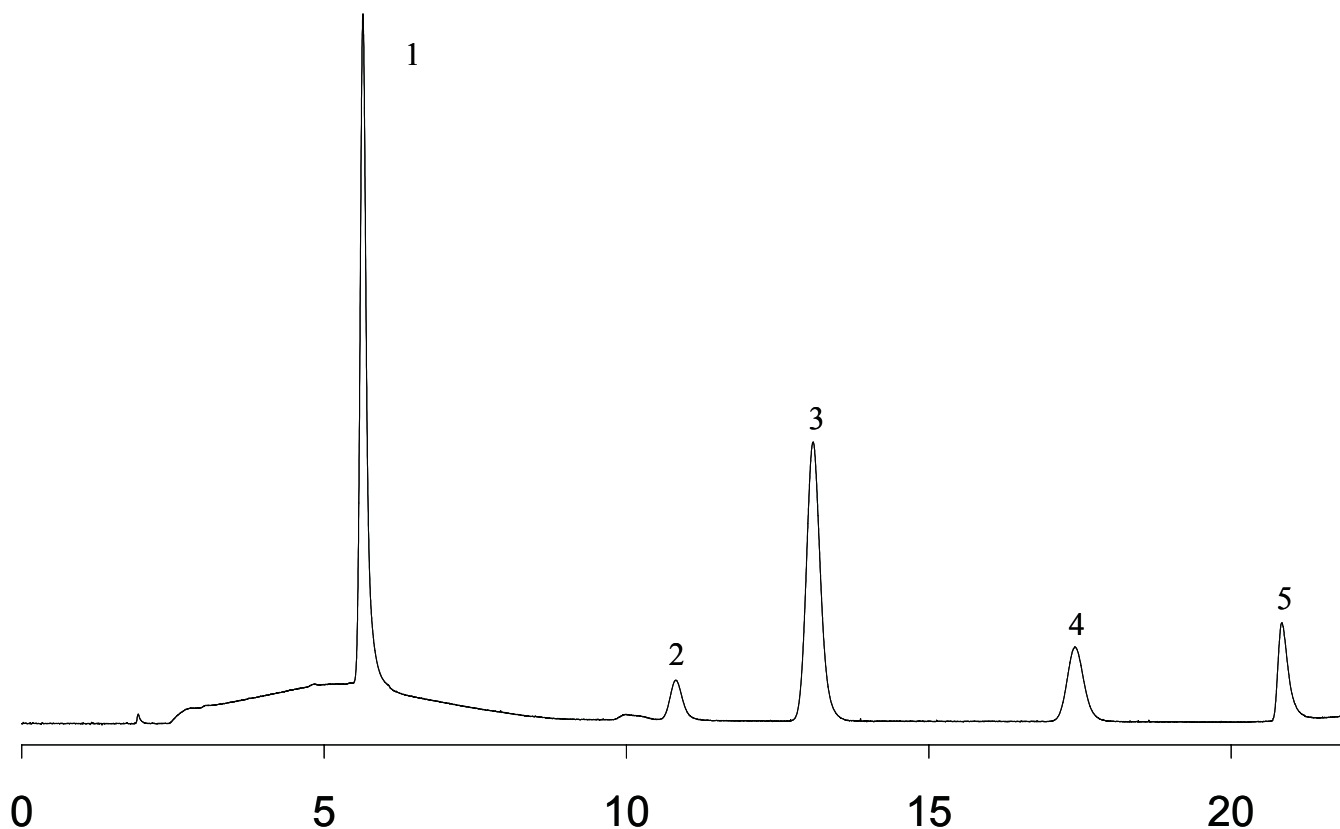

Phenomenex products are available worldwide. For the distributor in your country, contact Phenomenex USA, International Department by telephone, fax or e-mail: international@phenomenex.com.

 **phenomenex®**
 ...breaking with tradition™

USA	Puerto Rico	Canada	United Kingdom	Germany	New Zealand	Australia
mail: 411 Madrid Avenue Torrance, CA 90501-1430 USA	273 Sierra Morena, Suite #104 San Juan, Puerto Rico 00926	411 Madrid Avenue Torrance, CA 90501-1430 USA	Queens Avenue, Hurdfield Ind. Est., Macclesfield, Cheshire SK10 2BN, England	Zeppelinstr. 5 63741 Aschaffenburg Deutschland	P.O. Box 31-601 Milford Auckland New Zealand	P.O. Box 4084 Lane Cove, NSW 1595 Australia
tel.: (310) 212-0555	(800) 541-HPLC	(800) 543-3681	01625-501367	06021-58830-0	09-4780951	02-9428-6444
fax: (310) 328-7768	(310) 328-7768	(310) 328-7768	01625-501796	06021-58830-11	09-4780952	02-9428-6445
email: info@phenomenex.com	info@phenomenex.com	info@phenomenex.com	ukinfo@phenomenex.com	anfrage@phenomenex.com	info@phenomenex.co.nz	info@phenomenex.com.au