


Robitussin Maximum Strength on Synergi Hydro-RP

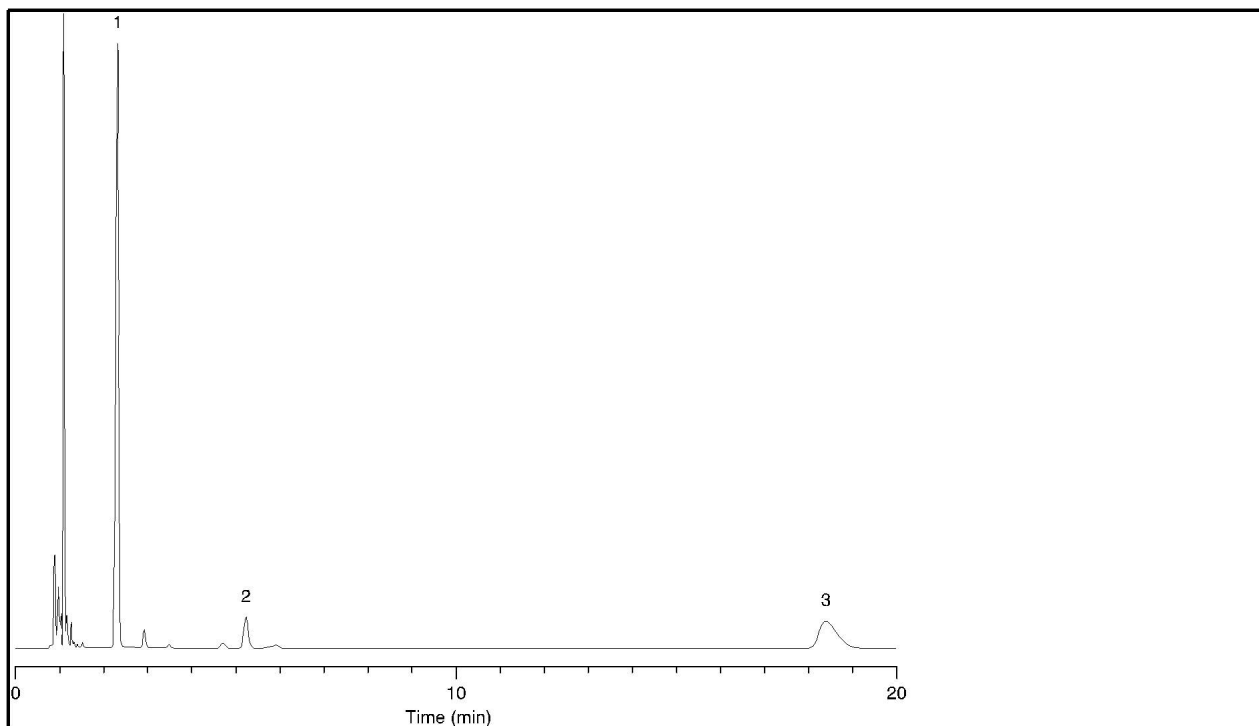

Column: Synergi 4 μ m C18 / ODS Hydro-RP

Dimensions: 150 x 4.6 mm ID

Order No:

Elution Type: Isocratic

Eluent A: 20mM KH₂PO₄ + 0.1% Hexane sulfonate, adjust to pH 3 **Pct. A:** 60

Eluent B: Acetonitrile **Pct. B:** 30

Eluent C: Methanol **Pct. C:** 10

Flow Rate: 1.5 mL/min

Temperature: 22 °C

Detection: UV-Vis Abs. - Variable Wave. @ 210 nm

SAMPLE ANALYTE

- 1 *Pseudoephedrine*
- 2 *Benzoic acid*
- 3 *Dextromethorphan*

SPE:

For more information call your Phenomenex Representative at (310) 212-0555

Copyright 2000 Phenomenex Inc.

Ref. No.: 14219