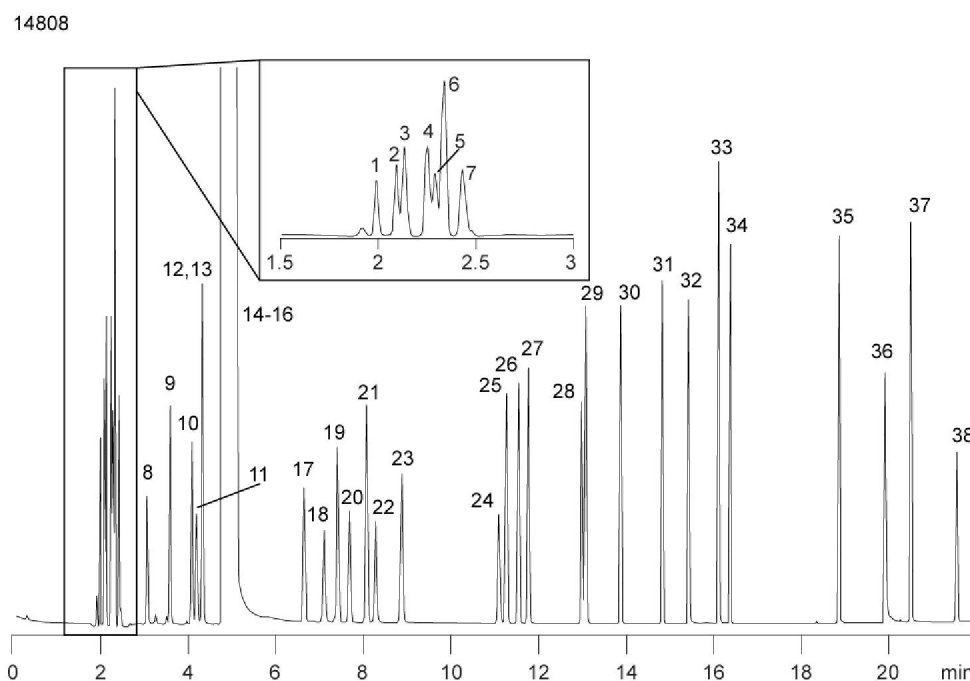


Industrial Solvents on ZB-WAX 30x0.25x0.25

Page 1 of 2

Column: Zebron ZB-WAX
Phase: Polyethylene glycol (PEG)
Dimensions: 30 meters x 0.25 mm x 0.25 µm
Order No: 7HG-G007-11
Oven Profile: 35°C for 5 min to 60°C at 4°C/min to 180°C at 12°C/min for 4 min.
Carrier Gas: Constant Flow Helium, 0.7 mL/min
Injection: split 100:1 1 µL @ 250 °C
Detection: Mass Selective (MSD) (180 °C)

Products used in this Application:

ANALYTES:

- | | | |
|---------------|-----------------------|------------------------|
| 1 Pentane | 10 Ethyl acetate | 19 Tetrachloroethylene |
| 2 Hexane | 11 Methanol | 20 Chloroform |
| 3 Ethyl ether | 12 t-Butanol | 21 Toluene |
| 4 Isooctane | 13 MEK | 22 1-Propanol |
| 5 MTBE | 14 Methylene chloride | 23 1,4-Dioxane |
| 6 Heptane | 15 Ethanol | 24 Allyl alcohol |
| 7 Cyclohexane | 16 Benzene | 25 Ethylbenzene |
| 8 Acetone | 17 Trichloroethylene | 26 p-Xylene |
| 9 THF | 18 Acetonitrile | 27 m-Xylene |

©Copyright 2000-2007 Phenomenex Inc.

For more information call your Phenomenex Representative at (310) 212-0555

www.phenomenex.com Phenomenex products are available worldwide. Email us at international@phenomenex.com					
Australia tel.: 02-9428-6444 email: info@phenomenex.com.au	Austria 01-319-1301 anfrage@phenomenex.com	Canada (800) 543-3681 info@phenomenex.com	Denmark 4824 8048 dkinfo@phenomenex.com	France 01 30 09 21 10 franceinfo@phenomenex.com	Germany 06021-58830-0 anfrage@phenomenex.com
Ireland tel.: 01 247 5405 email: eireinfo@phenomenex.com	Italy 051 736176 italiainfo@phenomenex.com	New Zealand 09-4780951 info@phenomenex.co.nz	Puerto Rico (800) 541-HPLC info@phenomenex.com	United Kingdom 01625-501367 ukinfo@phenomenex.com	USA (310) 212-0555 info@phenomenex.com

Industrial Solvents on ZB-WAX 30x0.25x0.25

Page 2 of 2

- 28 Pyridine
- 29 o-Xylene
- 30 Chlorobenzene
- 31 Styrene
- 32 3-Picoline
- 33 N,N-dimethylformamide
- 34 Anisole
- 35 Benzaldehyde
- 36 Ethylene glycol
- 37 N-methyl-2-Pyrrolidinone
- 38 Formamide

 www.phenomenex.com Phenomenex products are available worldwide. Email us at international@phenomenex.com					
Australia tel.: 02-9428-6444 email: info@phenomenex.com.au	Austria 01-319-1301 anfrage@phenomenex.com	Canada (800) 543-3681 info@phenomenex.com	Denmark 4824 8048 dkinfo@phenomenex.com	France 01 30 09 21 10 franceinfo@phenomenex.com	Germany 06021-58830-0 anfrage@phenomenex.com
Ireland tel.: 01 247 5405 email: eireinfo@phenomenex.com	Italy 051 736176 italiainfo@phenomenex.com	New Zealand 09-4780951 info@phenomenex.co.nz	Puerto Rico (800) 541-HPLC info@phenomenex.com	United Kingdom 01625-501367 ukinfo@phenomenex.com	USA (310) 212-0555 info@phenomenex.com