


Amoxicillin - Alternative Method


Column: Luna 5u C8(2) 100A
Dimensions: 150 x 4.6 mm ID
Order No: 00F-4249-E0
Elution Type: Isocratic
Eluent A: 20mM Ammonium acetate, pH 4.9
Eluent B: Methanol
Flow Rate: 1 mL/min
Col. Temp.: ambient
Detection: UV-Vis Abs. - Variable Wave. @ 230 nm (ambient)

Pct. A: 95**Pct. B:** 5

Page 1 of 1

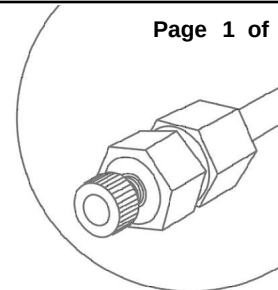

Products used in this Application:

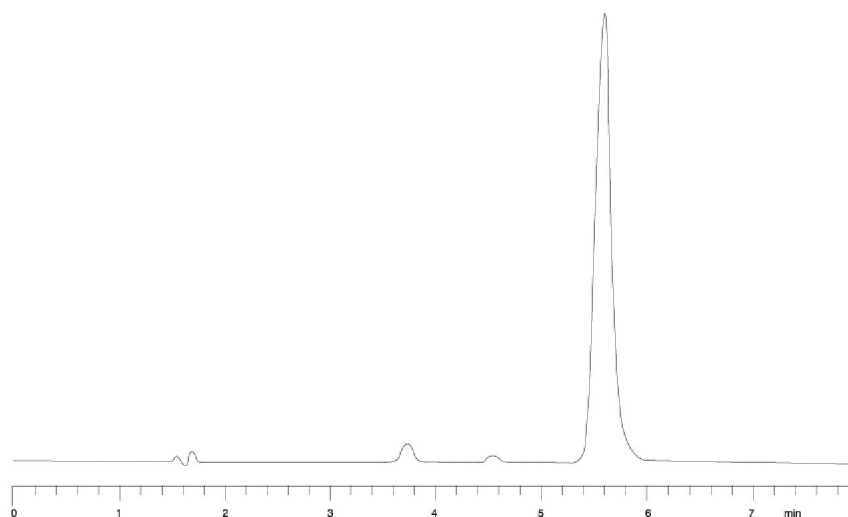LUNA[®]

ANALYTES:

1 Amoxicillin